

## Protocol

# Nuclear & Cytoplasmic Extract

The protocol for the isolation of nuclear and cytoplasmic fractions can be used to generate material for the gel shift/super shift/EMSA assays or any other assay requiring similar source material. Remember to use only specially prepared reagents that are free of any contaminating proteases or nucleases and to keep the preparation on ice where indicated. A common troubleshooting finding is that breakdown products caused by non-specific proteolysis are recognized in systems typically as lower molecular weight bands.

## Reagents Required

Product	Preparation	Item No.
Molecular Biology Grade Water		MB-010-0100
Cells for Nuclear Extract	Prepare cells from a cell line known to be positive for the intended assay. For NFkB HeLa or Jurkat cells can be used.	
Phosphate Buffered Saline (PBS)	Use 10X PBS, pH 7.2 (0.2 M Potassium Phosphate, 1.5 M NaCl). Dilute appropriate volume to 1X with molecular biology grade water.	MB-008
Cytoplasmic Extract (CE) Buffer with NP-40	Prepare a 1X solution composed of 10 mM HEPES, 60 mM KCl, 1 mM EDTA, 0.075% (v/v) NP-40, 1mM DTT and 1 mM PMSF, adjusted to pH 7.6. Convenient concentrated stocks of these reagents can be prepared such that 10 mL volumes of 1X CE buffer can be easily prepared.	
CE Buffer without Detergent	Similar to above buffer but without the addition of NP-40.	
Nuclear Extract (NE) Buffer	Prepare a 1X solution composed of 20 mM Tris Cl, 420 mM NaCl, 1.5 mM MgCl <sub>2</sub> , 0.2 mM EDTA, 1 mM PMSF and 25% (v/v) glycerol, adjusted to pH 8.0. Convenient concentrated stocks of these reagents can be prepared such that 10 mL volumes of 1X CE buffer can be easily prepared.	
5 M NaCl Solution	Prepare a solution of sodium chloride at 5.0 M.	
Glycerol		

## Procedure

1. Prepare approximately  $4 \times 10^7$  cells per extract.
2. Wash the cells gently with PBS buffer. Collect the cells by centrifugation using a microcentrifuge at 1000 rpms.
3. Resuspend the pellet in 5 pellet volumes of CE buffer (approximately 100  $\mu$ L).
4. Incubate on ice for 3 minutes to lyse the cells.

5. Spin the preparation using a microcentrifuge at 1000–1500 rpms for 4 minutes.
6. Move the supernatant containing the cytoplasmic extract to a clean tube (CE tube).
7. Wash the nuclei (NE tube) with 100  $\mu$ L of CE buffer without detergent.  
**Note:** Be careful to resuspend the fragile nuclei gently.
8. Spin the nuclei as above at 1000–1500 rpms for 4 minutes.
9. Remove the supernatant and add 1 pellet volume NE buffer to the nuclear pellet (approximately 50  $\mu$ L).
10. Adjust the salt concentration to 400 mM using 5 M NaCl (add ~35  $\mu$ L).
11. Add an additional pellet volume of NE buffer. Vortex to resuspend the pellet.
12. Incubate the extract on ice for 10 minutes. Vortex the mixture periodically to re-suspend the pellet.
13. Spin the CE and NE tubes at maximum speed for 10 minutes to pellet any nuclei.
14. Transfer the supernatant of the CE tube and NE tube separately to clean tubes. Add glycerol to the CE tube to 20%. Store at  $-70^{\circ}\text{C}$ .

---

Rockland products are for research use only and are not intended for therapeutic or diagnostic applications. Please contact a technical service representative for more information. All properties listed are typical characteristics and are not specifications. All suggestions and data are offered in good faith but without guarantee as conditions and methods of use of our products are beyond our control. All claims must be made within 30 days following the date of delivery. Suggested uses of our products are not recommendations to use our products in violation of any patent or as a license under any patent of Rockland Immunochemicals, Inc. If you require a commercial license to use this material and do not have one, then return this material, unopened to: Rockland Immunochemicals, Inc., P.O. BOX 5199, Limerick, Pennsylvania 19468, USA.

© 2025 Rockland Immunochemicals, Inc. All rights reserved.